

ANNEX I

METHOD FOR THE DETECTION AND DIAGNOSIS OF THE RING ROT BACTERIUM,
CLAVIBACTER MICHIGANENSIS (Smith) Davis et al. ssp. SEPEDONICUS
(Spieckermann et Kotthof) Davis et al. IN BATCHES OF POTATO TUBERS

Status: This is the original version (as it was originally adopted).

Appendix 6

MEDIA FOR GROWTH AND ISOLATION OF *C. SEPEDONICUM*

Yeast peptone glucose agar (YPGA)

Difco bacto yeast extract (No 0127)	5 g
Difco bacto peptone (No 0118)	5 g
D(+)-glucose (monohydrate)	10 g
Difco bacto purified agar (No 0560)	15 g
Distilled water	1 litre

Sterilize 0,5 litre volumes of medium by autoclaving at 115 °C for 20 minutes.