

## SCHEDULE 2

### TESTING METHODS

#### PART II

##### METHODS FOR THE ISOLATION OF *salmonella*

#### Day 4

29. The plates shall be examined and a minimum of 3 colonies from each plate showing suspicion of *Salmonella* growth shall be subcultured –

- (a) onto a nutrient agar plate;
- (b) onto a MacConkey agar plate<sup>(1)</sup>; and
- (c) into/onto biochemical media suitable for the identification of *Salmonella*.

These media shall be incubated at 37°C ± 1°C overnight.

---

(1) MacConkey aga – See (1963) International Standards for Drinking Water, World Health Organisation Distribution and Sales, CH-1211, Geneva 27, Switzerland