

## SCHEDULE 2

### Testing Methods

## PART 3

### METHOD FOR THE ISOLATION OF *ENTEROBACTERIACEAE*

#### **Samples with colonies of *Enterobacteriaceae***

6. After counting the colonies, characteristic colonies shall be taken at random from the agar plates, the number being at least the square root of the colonies counted. The colonies shall be subcultured onto a blood agar plate and incubated aerobically at  $37^{\circ}\text{C}\pm 1^{\circ}\text{C}$  for 20 hours $\pm$ 2 hours.